



Exciting News!

On July 15, 2019, SKC Inc. acquired Colormetric Laboratories, Inc. (CLI).

SKC is excited to add CLI surface and dermal assessment and control tools to our highly respected, long-standing sampling solutions that help you maintain the health and safety of workers. SKC looks forward to continuing to provide the same great CLI products with excellent customer and technical support. Learn more about SKC Inc. at <https://www.skcinc.com/about>.

Surface SWYPE for Acid/Base Cat. No. 769-1026

1. Lightly spray the area or item (e.g., workbench, tool, control knob) with Developing Solution for Acid/Base Cat. No. 769-1066.
2. Wait at least 30 seconds and then wipe with a Surface SWYPE pad.
3. The acid detector will change from neutral (orange) color to fuchsia/magenta at pH 3. For caustic exposures, the indicator will change to blue at pH 9.5.

SKC Limited Warranty and Return Policy

SKC products are subject to the SKC Limited Warranty and Return Policy, which provides SKC's sole liability and the buyer's exclusive remedy. To view the complete SKC Limited Warranty and Return Policy, go to <http://www.skcinc.com/warranty>.